

The logo for Mirus, featuring the word "Mirus" in a white serif font with a red checkmark above the 'i'. Below the text is a red wavy line.

**Mirus**

The Transfection Experts

A detailed 3D illustration of several cells. The cells are semi-transparent, revealing internal organelles like nuclei and mitochondria. One cell in the foreground is particularly large and shows a glowing orange nucleus with a DNA double helix structure. The background is dark with some greenish particles.


[www.TheTransfectionExperts.com](http://www.TheTransfectionExperts.com)

# THE TRANSFECTION EXPERTS

From basic research to product development, Mirus Bio scientists utilize their expertise in nucleic acid chemistry, cellular and molecular biology to establish innovative technologies and products for use in life science research. Over a decade ago, our first products reflected advancements achieved in our non-viral gene therapy research. Since then, our scientists have made significant contributions providing new and better tools for a variety of research applications in the life sciences.

Our technologies and expertise precipitated the acquisition of Mirus Bio Corporation's therapeutic division by Hoffman-La Roche in 2008. However, Mirus Bio's biological research products division remains as a stand-alone entity, called Mirus Bio LLC.

**At Mirus Bio LLC, we will continue to grow our expertise in nucleic acid delivery research and products to better serve you.**

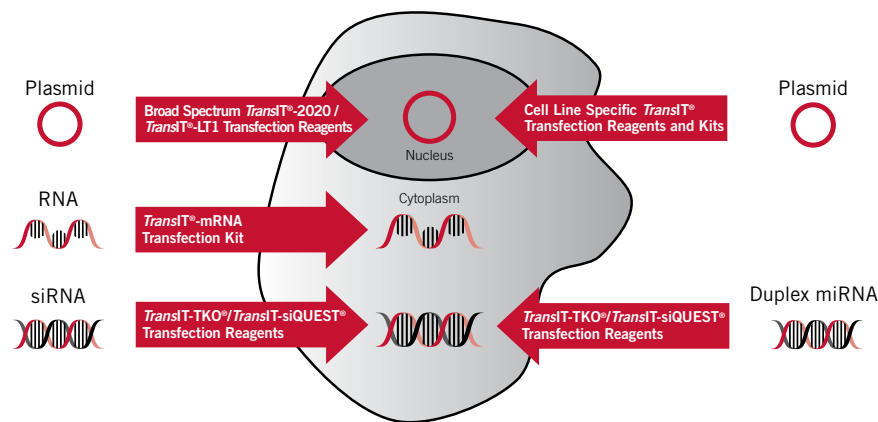

## WHAT IS TRANSFECTION?

Transfection is the introduction of any nucleic acid molecule by non-viral means into cultured mammalian cells. It is a common and convenient technology that aids in the study of gene function. Transfection methods can be divided into physical or direct transfer methods such as electroporation or chemical mediated transfection.

### Optimize Transfection Performance For All Reagents, Cells and Nucleic Acids

- 1. Ratio of transfection reagent to nucleic acid**  
In parallel for each cell type, transfect different amounts of nucleic acid with various amounts of transfection reagent. Titrate the amount of reagent while maintaining constant DNA concentration. Additionally, DNA concentration can be scaled.
- 2. Transfection complex formation time**  
Test different transfection complex formation times (reagent + nucleic acid incubation time) for each reagent and cell type being transfected.
- 3. Cell density (confluency) at time of transfection**  
Transfect each cell type at various densities ranging from 50% to 90% and monitor transfection performance.
- 4. Post-transfection incubation time (time after addition of transfection complexes to cells)**  
Harvest cells at different times post-transfection to determine the optimal assay time that meets your experimental goals.

For specific recommendations, please visit [www.mirusbio.com](http://www.mirusbio.com) or contact our transfection experts at [techsupport@mirusbio.com](mailto:techsupport@mirusbio.com).

# CONTENTS

## ▶ DNA TRANSFECTION

### ▷ BROAD SPECTRUM

<i>TransIT</i> ®-2020 Transfection Reagent.....	3
<i>TransIT</i> ®-LT1 Transfection Reagent .....	5

### ▷ CELL LINE SPECIFIC

<i>TransIT</i> ®-293 Transfection Reagent.....	7
<i>TransIT</i> ®-3T3 Transfection Kit .....	7
<i>TransIT</i> ®-CHO Transfection Kit.....	7
<i>TransIT</i> ®-COS Transfection Kit.....	7
<i>TransIT</i> -HeLaMONSTER® Transfection Kit.....	8
<i>TransIT</i> ®-Jurkat Transfection Reagent .....	8
<i>TransIT</i> ®-Keratinocyte Transfection Reagent .....	8
<i>TransIT</i> -Neural® Transfection Reagent.....	8
<i>TransIT</i> ®-Prostate Transfection Kit.....	8

### ▷ ELECTROPORATION

Ingenio™ Electroporation Kit .....	9
------------------------------------	---

### ▷ NEW! PROTEIN PRODUCTION

<i>TransIT</i> -PRO™ Transfection Kit .....	11
---	----

### ▷ PLASMID CONTROLS

Plasmid Delivery Control, Cy™3 .....	13
Plasmid Delivery Control, Fluorescein .....	13

## ▶ RNA TRANSFECTION

### ▷ siRNA

<i>TransIT</i> -TKO® Transfection Reagent.....	14-15
<i>TransIT</i> -siQUEST® Transfection Reagent .....	14-15
<i>TransIT</i> ® siPAK Kit .....	14-15
<i>TransIT</i> ® siPAK Plus Kit .....	14-15
Ingenio™ Electroporation Kits.....	16
Ingenio™ Electroporation Solution.....	16
Ingenio™ Electroporation Accessories.....	16
RNAi Delivery Control, Cy™3 .....	17
RNAi Delivery Control, Fluorescein .....	17

### ▷ LARGE RNA (VIRAL RNA and mRNA)

<i>TransIT</i> ®-mRNA Transfection Kit.....	18
---	----

## OUR PROMISE TO YOU: SERVICE THROUGH EXPERTISE

**Order Online 24 Hours a Day (Credit Cards or Purchase Orders)**

[www.mirusbio.com](http://www.mirusbio.com)

### Customer Service

Hours of Operation: 8:00 AM–4:00 PM Central Time, Monday–Friday

**email:** [sales@mirusbio.com](mailto:sales@mirusbio.com)

**phone:** 888.530.0801 (toll free within the U.S.) or +1.608.441.2852

**fax:** +1.608.441.2849

### Technical Service

Hours of Operation: 8:00 AM–4:00 PM Central Time, Monday–Friday

**email:** [techsupport@mirusbio.com](mailto:techsupport@mirusbio.com)

**phone:** 888.530.0801 (toll free within the U.S.) or +1.608.441.2852

**fax:** +1.608.441.2849

### Mail

Mirus Bio LLC

Attention: Customer Service

545 Science Drive

Madison, WI 53711 USA

### U.S. Distribution

Fisher Scientific



### International Distribution

Mirus Bio products are available worldwide through our international distributors. For a complete list, please visit:

[www.mirusbio.com/contact/distributors.aspx](http://www.mirusbio.com/contact/distributors.aspx) and contact the distributor in your region.

## Broad Spectrum

# TransIT®-2020 TRANSFECTION REAGENT

- ▷ **Superior Performance**—TransIT-2020 outperforms FuGENE® HD, Lipofectamine™ 2000, and Lipofectamine™ 2000 CD
- ▷ **Broad Spectrum DNA Delivery**—Achieve high expression in many cell types, including hard-to-transfect cells and insect cells
- ▷ **Animal Origin Free**—TransIT-2020 provides high performance with maximum compatibility


### PRODUCT SPECIFICATIONS

PRODUCT NO.	QUANTITY
MIR 5400	1.0 ml
MIR 5404	0.4 ml
MIR 5405	5 × 1.0 ml
MIR 5406	10 × 1.0 ml

To inquire about bulk pricing, please call 888.530.0801.

### Description


TransIT-2020 Reagent is a new and versatile transfection solution for broad spectrum DNA delivery into mammalian cells. This premium, broad spectrum reagent is animal component free allowing maximum compatibility for all downstream applications while outperforming major competitors in many cell types.

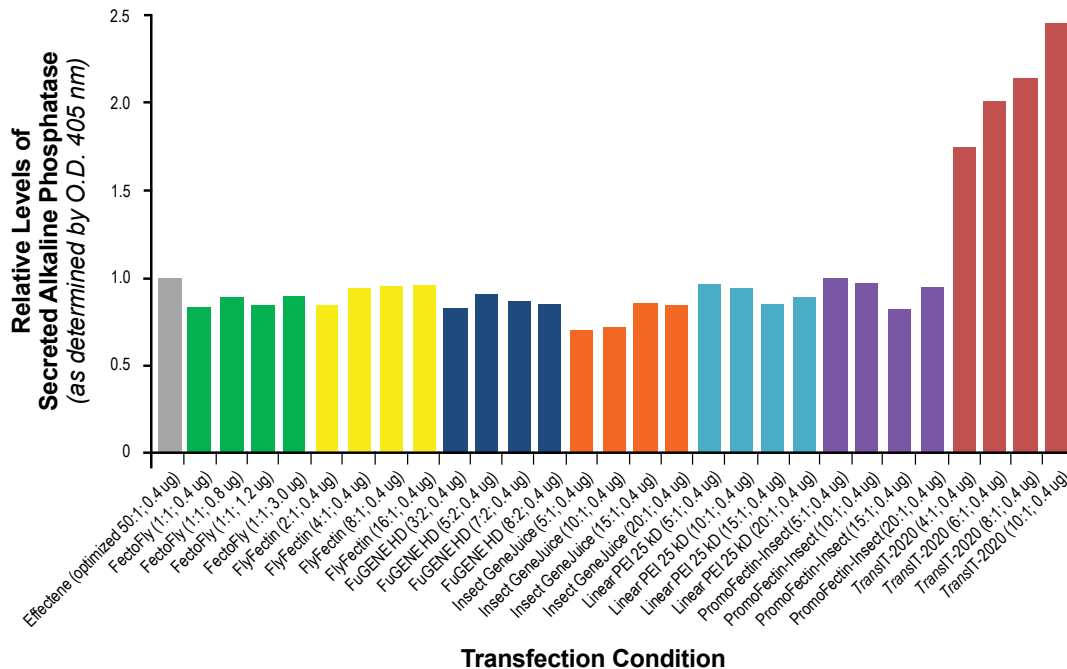

**FIGURE 1. The TransIT®-2020 Reagent Effectively Transfects *Drosophila* S2 Cells.** Cells were transfected with a plasmid construct expressing a secreted form of the Dscam extracellular domain fused to alkaline phosphatase (AP) in a 24-well plate. The ratio of transfection reagent to DNA and micrograms of DNA per well is noted beneath each bar. All products were used according to manufacturers' protocol. All transfections were performed in serum-free media for four hours followed by complete media supplementation. An AP enzymatic assay was used to measure the AP levels 24 hours post-transfection.

Data courtesy of Woj Wojtowicz, University of California, Berkeley

## **TransIT-2020®: High Performance, Animal free, Broad Spectrum DNA Transfection**

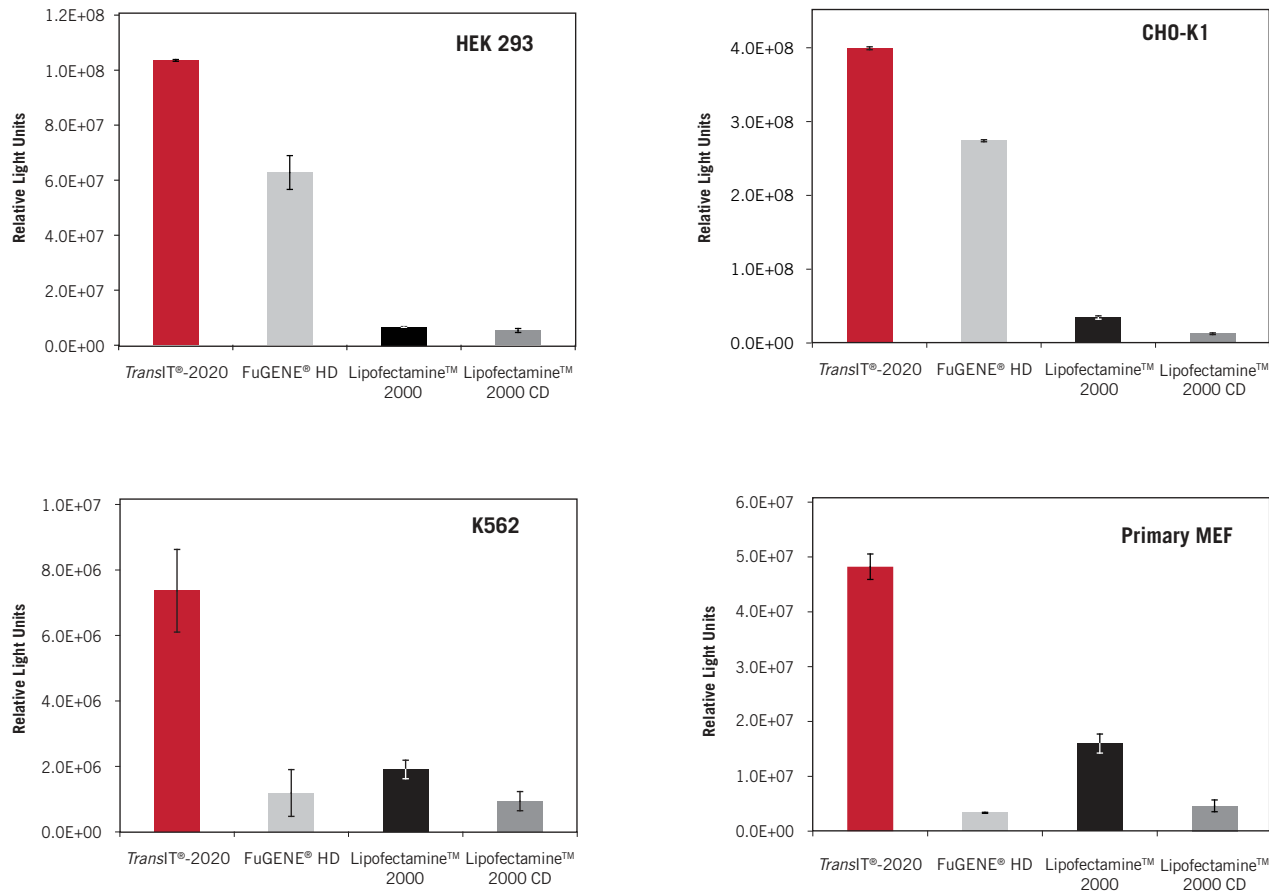

**FIGURE 2. Superior Gene Expression in a Broad Spectrum of Cell Types Using TransIT-2020® Transfection Reagent.** Luciferase gene expression was compared in HEK 293, K562, CHO-K1, and primary MEF cells transfected with a luciferase expression plasmid using TransIT®-2020 (Mirus Bio, 3:1 reagent to DNA ratio), FuGENE HD (Roche, 7:2 reagent to DNA ratio), Lipofectamine™ 2000 (Life Technologies Corporation, 5:2 reagent to DNA ratio), and Lipofectamine™ 2000 CD (Invitrogen, 5:2 reagent to DNA ratio). Transfections were performed in 24-well plates using 0.5 µg of plasmid DNA per well. The optimal level of each transfection reagent was determined empirically and all reagents were used according to manufacturer's protocol. Cells were harvested at 24 hours post-transfection and assayed for luciferase activity. Error bars represent the standard deviation of triplicate wells.

## Broad Spectrum

# TransIT®-LT1 TRANSFECTION REAGENT

- ▷ **Deliver Single or Multiple Plasmids**—Suitable for many applications such as siRNA/shRNA expression, gene expression, viral production, and promoter analysis
- ▷ **High Efficiency Delivery**—Achieve expression in a large population of cells to ensure experimental success
- ▷ **Low Cellular Toxicity**—Maintain cell density and reduce experimental biases due to toxicity induced cellular changes
- ▷ **Save Time**—No media changes or extensive optimization required
- ▷ **Save Money**—Utilize one transfection reagent and protocol for a variety of cells

## PRODUCT SPECIFICATIONS


PRODUCT NO.	QUANTITY
MIR 2304	0.4 ml
MIR 2300	1.0 ml
MIR 2305	5 × 1.0 ml
MIR 2306	10 × 1.0 ml

To inquire about bulk pricing, please call 888.530.0801.

### Description


TransIT®-LT1 Reagent is a broad spectrum, high efficiency DNA transfection reagent that is easy to use and exhibits minimal cellular toxicity. This reagent is a proprietary formulation of histone and cationic lipids that efficiently transfects cells in the presence or absence of serum.

**FIGURE 3. High Efficiency Transfection on a Broad Range of Cell Lines Using TransIT®-LT1 Reagent.** Various cell lines were transfected with an EGFP expression vector using TransIT®-LT1 Transfection Reagent in complete growth media with no media changes and analyzed by flow cytometry.

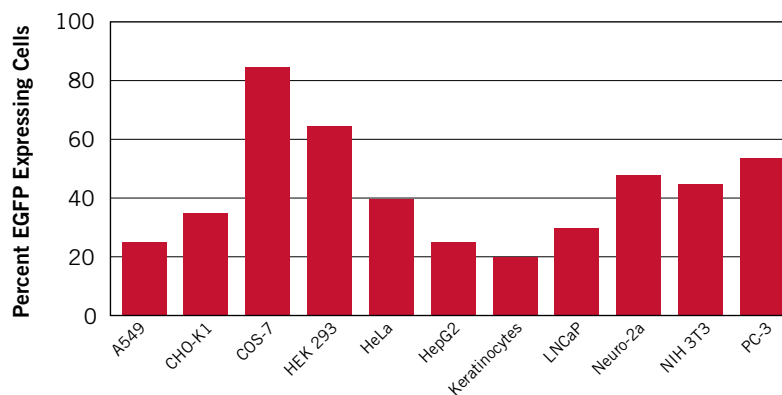

## *TransIT*<sup>®</sup>-LT1: High Efficiency, Low Toxicity Broad Spectrum Reagent

- ▷ **Physiologically relevant data you can trust**—Due to the exceptionally low cellular toxicity of *TransIT*-LT1 Reagent
- ▷ **Higher transfection efficiency than Lipofectamine<sup>™</sup> 2000**—At least twice the expression level regardless of cell density

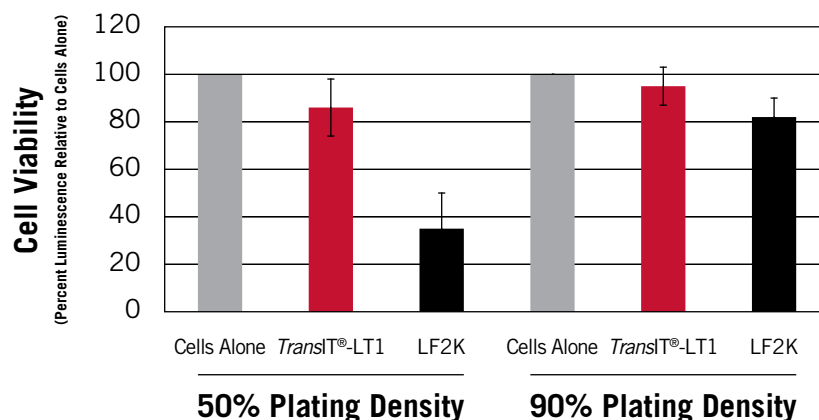

**FIGURE 4. High Cell Viability with *TransIT*<sup>®</sup>-LT1 Reagent, as Close to Nontransfected as Possible.** The impact of the *TransIT*<sup>®</sup>-LT1 and Lipofectamine<sup>™</sup> 2000 (LF2K) transfection reagents on cell viability was measured using the CellTiter-Glo<sup>®</sup> Luminescent Cell Viability Assay (Promega). This assay measures the levels of ATP in a cell population. HeLa cells at 50% and 90% confluency were transfected in duplicate using 3  $\mu$ l of *TransIT*<sup>®</sup>-LT1 Reagent or 4  $\mu$ l Lipofectamine<sup>™</sup> 2000 (LF2K) per well of a 12-well plate. Twenty-four hours post-transfection the cells were harvested and cell viability assayed. Data from three independent experiments performed on different days were averaged then scaled to untransfected HeLa controls (cells alone) and are represented as a percentage of the untransfected control.





## Cell Line Specific






# TransIT<sup>®</sup> CELL LINE SPECIFIC TRANSFECTION REAGENTS

The *TransIT*<sup>®</sup> Cell Line Specific DNA Transfection Reagents are formulated to maximize transfection efficiency while maintaining cellular health in many popular or hard-to-transfect cell lines types.

All of these reagents offer:

- ▷ **Unique Formulations**—Designed and optimized for each cell line or cell type to maximize transfection performance
- ▷ **Maximum Delivery Efficiency**—Enhanced transfection efficiency to ensure experimental success and improve experimental outcome
- ▷ **Low Cellular Toxicity**—Maintain cell density and reduce experimental biases due to toxicity induced cellular changes
- ▷ **Delivery of Single or Multiple Plasmids**—Suitable for many applications such as gene expression, siRNA/shRNA cassette expression, viral production and promoter analysis
- ▷ **Serum Compatible**—No media changes necessary or extensive optimization required, saving valuable research time

Product Name*	Applicable Cell Line(s) or Cell Type(s)	Efficiency**	Product No.	Quantity
<b>TransIT<sup>®</sup>-293 Transfection Reagent</b>				
	HEK 293, HEK 293T, and related	75–85%	MIR 2704	0.4 ml
			MIR 2700	1.0 ml
			MIR 2705	5 × 1.0 ml
			MIR 2706	10 × 1.0 ml
<b>TransIT<sup>®</sup>-3T3 Transfection Kit (TransIT<sup>®</sup>-3T3 Reagent and 3T3 Authority Reagent)</b>				
	NIH 3T3, 3T3-L1 and related	55–65%	MIR 2184	0.4 ml
			MIR 2180	1.0 ml
			MIR 2185	5 × 1.0 ml
			MIR 2186	10 × 1.0 ml
<b>TransIT<sup>®</sup>-CHO Transfection Kit (TransIT<sup>®</sup>-CHO Reagent and CHO Mojo Reagent)</b>				
	CHO-K1 and related	50–60%	MIR 2174	0.4 ml
			MIR 2170	1.0 ml
			MIR 2175	5 × 1.0 ml
			MIR 2176	10 × 1.0 ml
<b>TransIT<sup>®</sup>-COS Transfection Kit (TransIT<sup>®</sup>-COS Reagent and COS Boss Reagent)</b>				
	COS-1, COS-7 and related	90–95%	MIR 2194	0.4 ml
			MIR 2190	1.0 ml
			MIR 2195	5 × 1.0 ml
			MIR 2196	10 × 1.0 ml

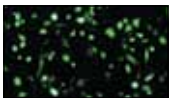Product Name*	Applicable Cell Line(s) or Cell Type(s)	Efficiency**	Product No.	Quantity
<b>TransIT<sup>®</sup>-HeLaMONSTER<sup>®</sup> Transfection Kit (TransIT<sup>®</sup>-HeLa Reagent and MONSTER Reagent)</b>				
	HeLa and related	50–60%	MIR 2904	0.4 ml
			MIR 2900	1.0 ml
			MIR 2905	5 × 1.0 ml
			MIR 2906	10 × 1.0 ml
<b>TransIT<sup>®</sup>-Jurkat Transfection Reagent</b>				
	Jurkat, Jurkat-E6, RAW 264.7, THP-1, K562, and other lymphoid cell lines	10–25%	MIR 2124	0.4 ml
			MIR 2120	1.0 ml
			MIR 2125	5 × 1.0 ml
			MIR 2126	10 × 1.0 ml
<b>TransIT<sup>®</sup>-Keratinocyte Transfection Reagent</b>				
	NIKS (Near Diploid Immortalized Keratinocytes) and primary keratinocytes	20–30%	MIR 2804	0.4 ml
			MIR 2800	1.0 ml
			MIR 2805	5 × 1.0 ml
			MIR 2806	10 × 1.0 ml
<b>TransIT<sup>®</sup>-Neural<sup>®</sup> Transfection Reagent</b>				
	C6, Daoy, DB-TRG-05MG, DI-TNC1, HCN-1A, Neuro-2a, PC-12, SK-N-MC, SVGp12	≥75%	MIR 2144	0.4 ml
			MIR 2140	1.0 ml
			MIR 2145	5 × 1.0 ml
			MIR 2146	10 × 1.0 ml
<b>TransIT<sup>®</sup>-Prostate Transfection Kit (TransIT<sup>®</sup>-Prostate Reagent and Prostate Boost Reagent)</b>				
	DU 145, LNCAP, PC3	>50%	MIR 2134	0.4 ml
			MIR 2130	1.0 ml
			MIR 2135	5 × 1.0 ml
			MIR 2136	10 × 1.0 ml

\* Single tube reagents contain the indicated transfection reagent. Transfection reagents with two components are named "Kits" and both components are listed following the product name.

\*\* Transfection efficiency determined by transfection of an EGFP expression vector followed by quantification of the percentage of cells expressing EGFP.

To inquire about bulk pricing, please call 888.530.0801.

## Electroporation

# INGENIO™ ELECTROPORATION PRODUCTS

- ▷ **High Efficiency Electroporation of Hard-to-Transfect Cells**—Conduct research in biologically relevant cells
- ▷ **Compatible with All Electroporation Instruments**—Use your existing system including an amaxa Nucleofector® or Bio-Rad Gene Pulser; no need to purchase additional specialized equipment
- ▷ **Save Money**—Replace your amaxa Nucleofector® Kits with the Ingenio™ Electroporation Kit and realize significant savings without sacrificing performance
- ▷ **Buy Only What You Need**—Ingenio Electroporation Solution is available alone or as part of a complete kit with cuvettes and cell droppers
- ▷ **Higher Cell Viability**—Less harmful than other electroporation methods

### Description

The Ingenio™ Electroporation Solution facilitates efficient and reliable delivery of nucleic acids to eukaryotic cells traditionally resistant to chemical transfection. Ingenio™ is a broad spectrum solution that supports high efficiency electroporation with minimal toxicity. It replaces standard electroporation solutions including phosphate buffered saline and serum-free media. Ingenio™ is compatible with multiple instruments and facilitates a wide range of applications requiring nucleic acid delivery to cells. The Ingenio Solution is available alone and as part of a kit with cuvettes and cell droppers.

### PRODUCT SPECIFICATIONS

#### Ingenio™ Electroporation Kits (solution, 0.4 cm cuvettes, cell droppers)

PRODUCT NO.	SIZE
MIR 50113	25 RXN <sup>a</sup>
MIR 50116	50 RXN <sup>a</sup>
MIR 50119	100 RXN <sup>a</sup>

#### Ingenio™ Electroporation Kits (solution, 0.2 cm cuvettes, cell droppers)

PRODUCT NO.	SIZE
MIR 50112	25 RXN <sup>a</sup>
MIR 50115	50 RXN <sup>a</sup>
MIR 50118	100 RXN <sup>a</sup>

#### Ingenio™ Electroporation Solution

PRODUCT NO.	SIZE	QUANTITY
MIR 50111	25 RXN <sup>b</sup>	6.25 ml
MIR 50114	50 RXN <sup>b</sup>	12.5 ml
MIR 50117	100 RXN <sup>b</sup>	25 ml


#### Ingenio™ Electroporation Accessories

PRODUCT NO.	DESCRIPTION	SIZE
MIR 50120	0.2 cm Cuvettes	25 PK
MIR 50121	0.2 cm Cuvettes	50 PK
MIR 50122	0.4 cm Cuvettes	25 PK
MIR 50123	0.4 cm Cuvettes	50 PK
MIR 50124	Cell Droppers	25 PK
MIR 50125	Cell Droppers	50 PK

<sup>a</sup> Electroporations per kit.


<sup>b</sup> Number of electroporations in 0.4 cm cuvette.

To inquire about bulk pricing, please call 888.530.0801.

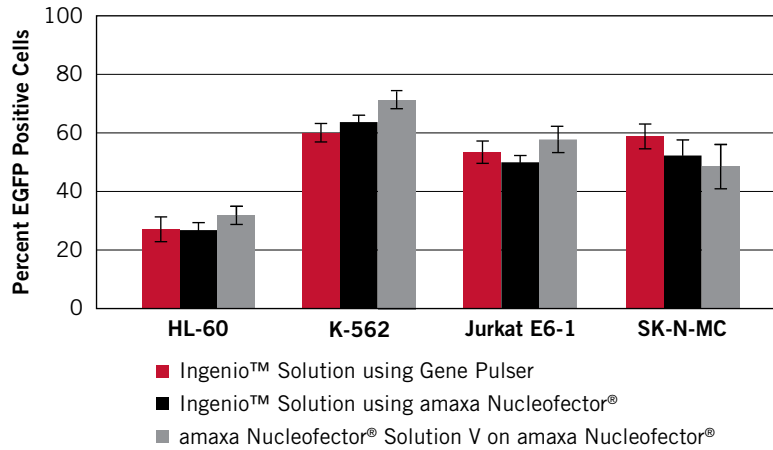

**FIGURE 5.** Ingenio™ Solution Provides Comparable Efficiency on amaxa's Nucleofector® Device. Cells were electroporated in parallel with an EGFP reporter vector. Two electroporators were used with different electroporation kits: the Ingenio™ Electroporation Kit was used in the Gene Pulser Xcell™ Eukaryotic System (Bio-Rad) and the amaxa Nucleofector™ II Device (Lonza); the amaxa Nucleofector Kit V was used in the amaxa Nucleofector™ II Device, all according to manufacturer's recommendations. EGFP expressing cells were identified 24 hours post-electroporation by flow cytometry and presented as a percentage of the live cell population. Experiments were performed in triplicate on three separate days and the data averaged.

**FIGURE 6.** Ingenio™ Electroporation Kits were used to transfect representative cell types using amaxa Nucleofector® Device. Cells were assayed at 24 hours post-electroporation by flow cytometry and reported as percentage of live cell population. Visit [www.TheTransfectionExperts.com](http://www.TheTransfectionExperts.com) for ideal pulse conditions.

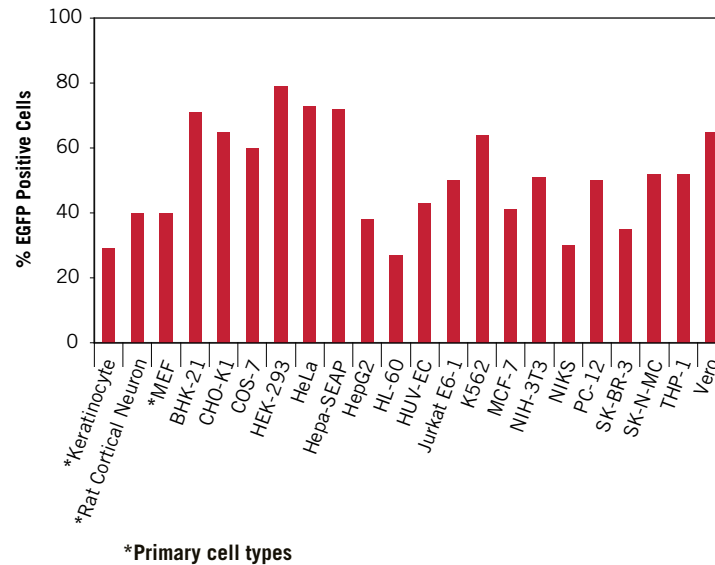

## Protein Production

# TransIT-PRO™ TRANSFECTION KIT

- ▷ **High Performance**—A cost-effective alternative for protein production
- ▷ **Easy to Use**—Compatible with multiple CHO media formulations
- ▷ **Reliable Scale-up**—Linear from 4 to 400 milliliters

### Description

Decrease time to produce usable protein by maximizing target protein yields through transient transfection. The *TransIT-PRO*™ Transfection Kit uses animal origin free components designed for high and reproducible protein yield in suspension CHO and 293 derived cells. Since it is compatible with varied media formulations, the same media can be used for both transient and stable expression. The *TransIT-PRO*™ outperforms linear PEI in protein yield, while providing a cost-effective alternative to FreeStyle™ MAX.


### PRODUCT SPECIFICATIONS

#### *TransIT-PRO*™ Transfection Kit

PRODUCT NO.	SIZE
MIR 5700	1 ml
MIR 5760	10 ml


To inquire about bulk pricing, please call 888.530.0801.

## High Performance: A Cost-effective Alternative for Protein Production

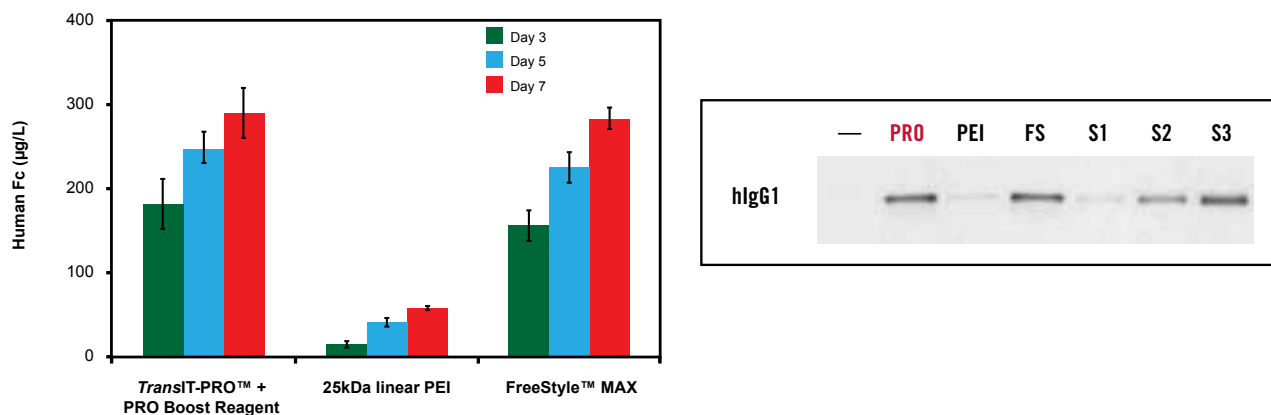


**FIGURE 7. Achieve High Antibody Titers Using the *TransIT-PRO*™ Transfection Kit.** Human IgG1 was produced by transient transfection using *TransIT-PRO* and PRO Boost Reagent (1:1:1), 25kDa linear PEI (6:1) or FreeStyle™ MAX (1:1) transfection reagents according to the manufacturers or published protocol (reagent:DNA ratio). Transfections were performed using 1 µg plasmid DNA per milliliter of culture and  $0.5 \times 10^6$  cells/ml at the time of transfection. FreeStyle™ CHO-S cells were cultured in 20ml of FreeStyle™ CHO Expression medium in 125 ml shake flasks. (A) Day 3, 5 and 7 supernatants were clarified and analyzed using a human IgG-Fc sandwich ELISA. Error bars represent the standard deviation of triplicate technical replicates, 25kDa linear PEI is duplicate technical replicates. (B) Day 7 supernatants were clarified and analyzed by Western blot. An IgG standard was included for quantification estimate (S1= 1.6 mg/L, S2= 3.2 mg/L, S3 = 6.3 mg/L).

## Easy to Use: Compatible with Multiple Media Formulations

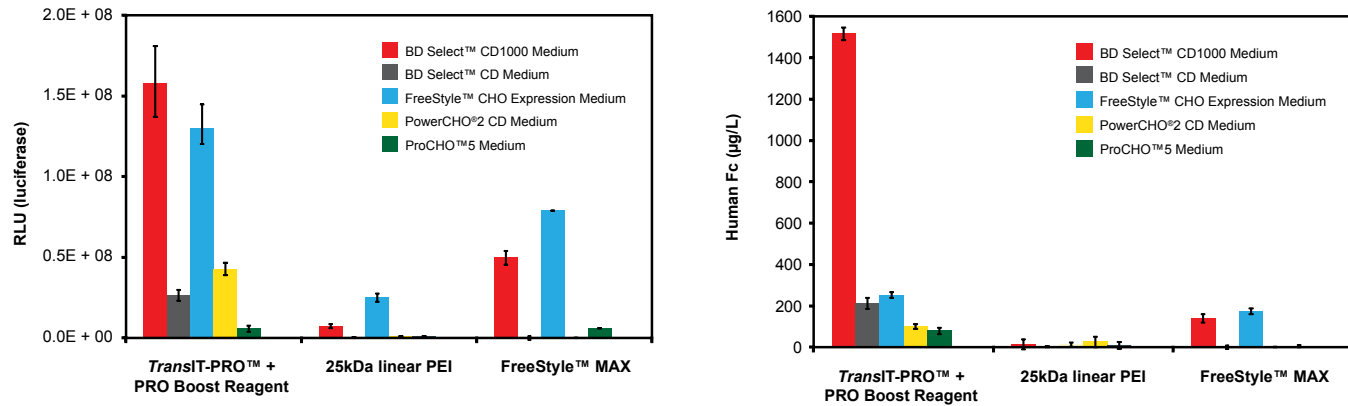

**FIGURE 8.** *TransIT-PRO™ Provides High Performance Across Varied Media Formulations.* FreeStyle™ CHO-S cells were adapted to five representative growth media including: BD Select™ CD1000 medium (Becton, Dickinson and Company, Franklin Lakes, NJ), BD Select™ CHO medium (Becton, Dickinson and Company, Franklin Lakes, NJ), FreeStyle™ CHO Expression medium (Life Technologies Corporation, Carlsbad, CA), ProCHO™5 medium (Lonza Inc., Allendale, NJ) and PowerCHO®2 CD medium (Lonza Inc., Allendale, NJ). Cells were transfected with a plasmid using the *TransIT-PRO* and PRO Boost Reagent (1:1:1), 25kDa linear PEI (4:1) (Polysciences, Warrington, PA), or FreeStyle™ Max (1:1) (Life Technologies Corporation, Carlsbad, CA) transfection reagents according to the manufacturers or published protocol (reagent:DNA ratio). Transfections were performed in 24-well deep well shaker blocks using 1 µg plasmid DNA per milliliter of culture and  $0.5 \times 10^6$  cells/ml at the time of transfection, 2 ml total culture volume. (A) Luciferase expression was assessed 48 hours post-transfection using a conventional luciferase assay. Error bars represent the standard deviation of duplicate wells. (B) Human IgG1 was quantitated from day 5 clarified supernatants and analyzed by a human anti-Fc sandwich ELISA. Error bars represent the standard deviation of triplicate wells.

## Scale-up: No Problem

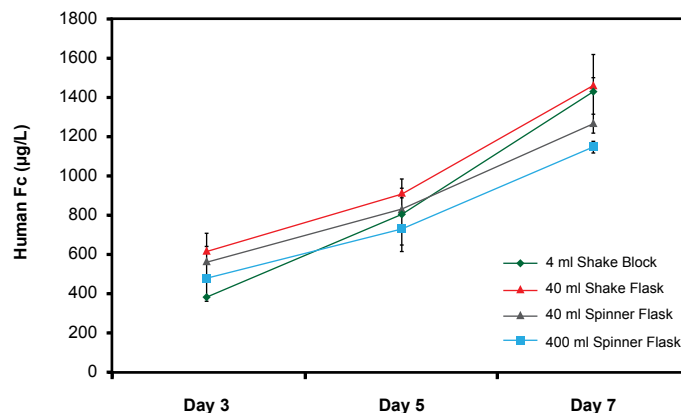

**FIGURE 9.** *Scaling of Transient Transfection Using TransIT-PRO™ Transfection Kit is Linear From 4 to 400 Milliliters.* Human IgG1 was produced by transient transfection using the *TransIT-PRO™* Transfection Kit and a plasmid encoding a human IgG1 construct. A DNA concentration was 1 µg/ml of culture and a ratio of *TransIT-PRO*:PRO Boost Reagent:DNA ratio of 1:0.5:1. Cells were plated at a density of  $0.5 \times 10^6$  cells/ml at the time of transfection. CHO-S cells were cultured in BD Select™ CD1000 media using 4 ml per well of a 24-well deep well shake block, 40 ml in 125 ml Erlenmeyer shake flask, 40 ml in 125 ml 2 sidearm spinner flask and 400 ml in 500 ml 2 sidearm spinner flask. Day 3, 5 and 7 supernatants were clarified and analyzed by an anti-Fc sandwich ELISA. Error bars represent the standard deviation of triplicate technical replicates.

## Plasmid Controls

### Label IT<sup>®</sup> PLASMID DELIVERY CONTROLS

- ▷ **Sensitive**—Easily visualize transfected cells and assess delivery efficiency using fluorescent microscopy
- ▷ **Compatible**—Suitable for co-delivery experiments with functional plasmids
- ▷ **Easy to Use**—Supplied as a ready to use 0.5 mg/ml solution for both *in vitro* and *in vivo* tracking studies

#### Description


The *Label IT*<sup>®</sup> Plasmid Delivery Controls consist of either Cy<sup>™</sup>3 or fluorescein labeled 2.7 kb plasmid for assessment of delivery efficiency in mammalian cells.

#### PRODUCT SPECIFICATIONS


##### *Label IT*<sup>®</sup> Plasmid Delivery Controls

LABEL	PRODUCT NO.	QUANTITY
Cy <sup>™</sup> 3	MIR 7904	10 µg
	MIR 7905	100 µg
Fluorescein	MIR 7906	10 µg
	MIR 7907	100 µg

To inquire about bulk pricing, please call 888.530.0801.

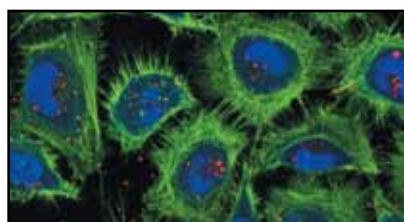

**FIGURE 10.** The *Label IT*<sup>®</sup> Cy<sup>™</sup>3 Plasmid Delivery Control Allows Quick Assessment of Delivery Efficiency. HeLa cells were transfected in complete media with the *Label IT*<sup>®</sup> Cy<sup>™</sup>3 Plasmid Delivery Control (red) using the *TransIT*<sup>®</sup>-LT1 Transfection Reagent. Twenty-four hours post-transfection, the cells were fixed, then counterstained to locate the nuclei (blue) and actin (green).

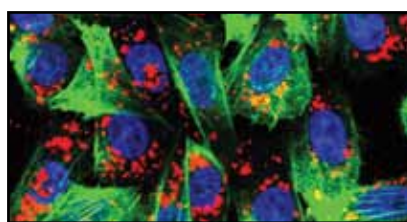

**FIGURE 11.** *Label IT*<sup>®</sup> Cy<sup>™</sup>3 Plasmid Delivery Control Allows Quick Assessment of Electroporation Efficiency. CHO-K1 cells were electroporated in Ingenio Electroporation Solution with *Label IT*<sup>®</sup> Cy<sup>™</sup>3 Plasmid Delivery Control (red). Twenty-four hours post-electroporation, cells were washed, fixed and counterstained to identify the nuclei (blue) and actin (green). The image was acquired using a confocal microscope.


## siRNA

# TransIT-TKO® & TransIT-siQUEST® TRANSFECTION REAGENTS

- ▷ **Broad Spectrum siRNA and Duplex miRNA Delivery**—Utilize each transfection reagent and protocol for a variety of cells
- ▷ **Two Different Reagent Formulations**—Two choices when identifying the best transfection reagent to maximize gene knockdown in a given cell line
- ▷ **High Knockdown Efficiency**—Achieve optimal gene silencing in a large percentage of cells to ensure experimental success
- ▷ **Low Cellular Toxicity**—Maintain cell density and reduce experimental biases due to alterations in cellular health
- ▷ **Flexible Protocol**—use with either standard or reverse transfections

### Description

*TransIT-TKO*® and *TransIT-siQUEST*® small RNA (siRNA and duplex miRNA) Transfection Reagents are broad spectrum reagents that are easy to use and exhibit minimal cellular toxicity. Each reagent is uniquely formulated and exhibits distinct siRNA/miRNA transfection profiles. These two reagents allow the user to identify the best siRNA and duplex

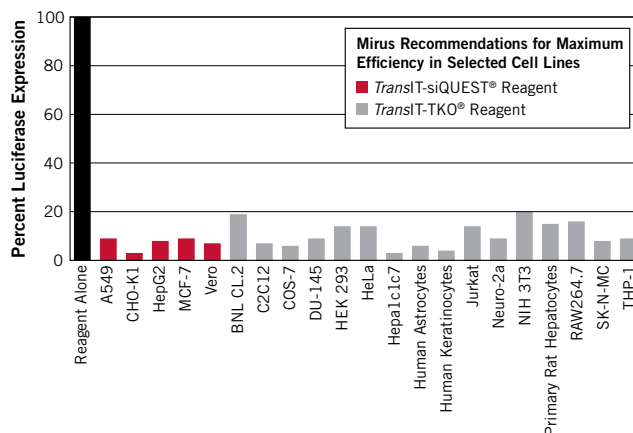

**FIGURE 12. Highly Efficient Knockdown of Transiently Expressed Firefly Luciferase.** Firefly and sea pansy luciferase reporter vectors were co-transfected into various cell lines. Subsequently, firefly luciferase expression was knocked down by transfection of 25 nM anti-firefly luciferase siRNA using either the *TransIT-siQUEST*® (red) or *TransIT-TKO*® (gray) Transfection Reagent. Bars indicate the percent of normalized firefly luciferase expression as compared to the reagent alone control 24 hours post-transfection.

miRNA transfection reagent for their particular cell line.

### PRODUCT SPECIFICATIONS

#### *TransIT-siQUEST*® Transfection Reagent

PRODUCT NO.	QUANTITY
MIR 2114	0.4 ml
MIR 2110	1.0 ml
MIR 2115	5 × 1.0 ml
MIR 2116	10 × 1.0 ml

#### *TransIT-TKO*® Transfection Reagent

PRODUCT NO.	QUANTITY
MIR 2154	0.4 ml
MIR 2150	1.0 ml
MIR 2155	5 × 1.0 ml
MIR 2156	10 × 1.0 ml

#### *TransIT*®-siPAK Kit

Trial sizes of *TransIT-TKO*® and *TransIT-siQUEST*® Transfection Reagents.

PRODUCT NO.	QUANTITY
MIR 2260	0.1 ml Each


#### *TransIT*®-siPAK Plus Kit

The *TransIT*®-siPAK with the addition of 10 µg the *Label IT*® RNAi Delivery Control, Fluorescein (MIR 7902).

PRODUCT NO.	QUANTITY
MIR 2270	0.1 ml Each

To inquire about bulk pricing, please call 888.530.0801.

## TransIT-TKO® & TransIT-siQUEST® continued

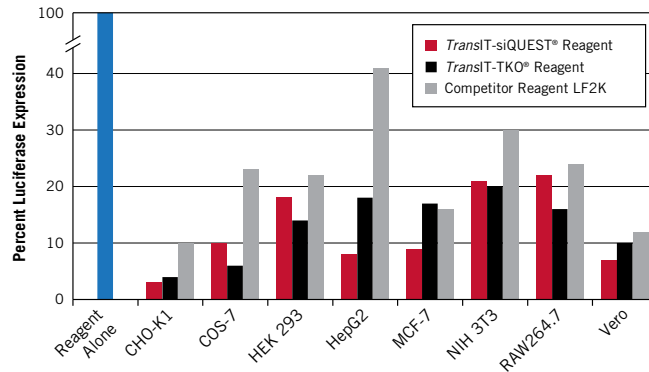

**FIGURE 13.** Knockdown Efficiencies Using *TransIT-siQUEST®* and *TransIT-TKO®* Reagents and Lipofectamine™ 2000 (LF2K). Firefly and sea pansy luciferase reporter vectors were co-transfected into various cell lines using the *TransIT®*-LT1 Reagent. Subsequently, firefly luciferase expression was knocked down by transfection of 25 nM anti-firefly luciferase siRNA using either *TransIT-siQUEST®* (red), *TransIT-TKO®* (black) or LF2K (gray) Reagents. Bars indicate the percent of normalized firefly luciferase expression as compared to each reagent alone control 24 hours post-transfection.

Cell Line (Source)	Endogenous Transcript	<i>TransIT-TKO®</i> Knockdown Efficiency	<i>TransIT-siQUEST®</i> Knockdown Efficiency
A549-luc (human lung)	Luciferase*	77%	82%
BNL CL.2 (mouse liver)	MAPK1 MAPK3	80% 83%	-- --
CHO-luc (hamster ovary)	Luciferase*	86%	91%
HEK 293-luc (human kidney)	Luciferase*	83%	77%
HeLa (human cervix)	Lamin A/C GAPDH	80% 80%	-- --
HeLa-luc (human cervix)	Luciferase*	84%	82%
Hepa-luc (mouse liver)	Luciferase*	--	92%
HepG2 (human liver)	MAPK1	80%	--
NIH 3T3-luc (mouse fibroblast)	Luciferase*	85%	89%
NIH 3T3-L1	MAPK1 MAPK3	70% 70%	-- --
Secondary Human Astrocytes	Lamin A/C	80%	--
Primary Mouse Hepatocytes	ABC A1 Lamin A/C PPAR-alpha	70% 81% --	-- -- 82%

**TABLE 1.** Knockdown of Endogenous Genes Using *TransIT-TKO®* or *TransIT-siQUEST®* Transfection Reagents. Cells were transfected with siRNAs targeting the indicated genes using the *TransIT-TKO®* or *TransIT-siQUEST®* Reagents, and the knockdown percentage was determined using quantitative RT-PCR or luciferase assays.

\* Firefly luciferase expression vectors were stably integrated into the parent cell lines and clonal lines constitutively expressing firefly luciferase were used.


## siRNA

# INGENIO™ ELECTROPORATION PRODUCTS

- ▷ **High Efficiency Electroporation of Hard to Transfect Cells**—Conduct research in biologically relevant cells
- ▷ **Compatible with All Electroporation Instruments**—Use your existing system including an amaxa Nucleofector® or Bio-Rad Gene Pulser; no need to purchase additional specialized equipment
- ▷ **Save Money**—Replace your amaxa Nucleofector® Kits with the Ingenio™ Electroporation Kit and realize significant savings without sacrificing performance
- ▷ **Buy Only What You Need**—Ingenio Electroporation Solution is available alone or as part of a complete kit with cuvettes and cell droppers
- ▷ **Higher Cell Viability**—Less harmful than other electroporation methods

### Description

The Ingenio™ Electroporation Solution facilitates efficient and reliable delivery of nucleic acids to eukaryotic cells traditionally resistant to chemical transfection. Ingenio™ is a broad spectrum solution that supports high efficiency electroporation with minimal toxicity. It replaces standard electroporation solutions including phosphate buffered saline and serum-free media. Ingenio™ is compatible with multiple instruments and facilitates a wide range of applications requiring nucleic acid delivery to cells. The Ingenio Solution is available alone and as part of a kit with cuvettes and cell droppers.

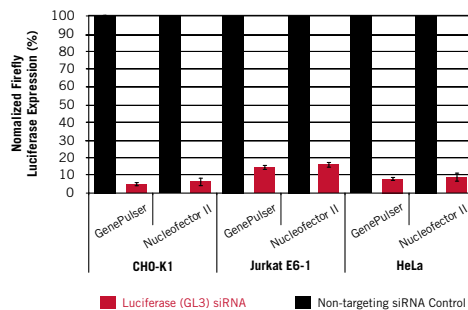

**FIGURE 14. The Ingenio™ Kits Effectively Electroporate siRNA on the amaxa Nucleofector® and the Gene Pulser® Electroporators.** siRNA and plasmid DNA were co-electroporated with the indicated cell lines with the Ingenio™ electroporation solution in 0.2 cm cuvettes using either the Gene Pulser® (Bio-Rad) or the amaxa Nucleofector® II (Lonza). 10µg/ml of plasmid encoding Firefly luciferase was co-electroporated with 250nM of either non-targeting siRNA control (Dharmacon Non-Targeting siRNA #1) or GL3 siRNA into Jurkat E6-1, HeLa and CHO-K1 cells. Twenty-four hours post electroporation, cells were harvested and assayed for luciferase activity. Data from independent experiments performed on different days were averaged then scaled to non-targeting siRNA control and are represented as a percentage of the control.

### PRODUCT SPECIFICATIONS

#### Ingenio™ Electroporation Kits (solution, 0.4 cm cuvettes, cell droppers)

PRODUCT NO.	SIZE
MIR 50113	25 RXN <sup>a</sup>
MIR 50116	50 RXN <sup>a</sup>
MIR 50119	100 RXN <sup>a</sup>

#### Ingenio™ Electroporation Kits (solution, 0.2 cm cuvettes, cell droppers)

PRODUCT NO.	SIZE
MIR 50112	25 RXN <sup>a</sup>
MIR 50115	50 RXN <sup>a</sup>
MIR 50118	100 RXN <sup>a</sup>

#### Ingenio™ Electroporation Solution

PRODUCT NO.	SIZE	QUANTITY
MIR 50111	25 RXN <sup>b</sup>	6.25 ml
MIR 50114	50 RXN <sup>b</sup>	12.5 ml
MIR 50117	100 RXN <sup>b</sup>	25 ml

#### Ingenio™ Electroporation Accessories

PRODUCT NO.	DESCRIPTION	SIZE
MIR 50120	0.2 cm Cuvettes	25 PK
MIR 50121	0.2 cm Cuvettes	50 PK
MIR 50122	0.4 cm Cuvettes	25 PK
MIR 50123	0.4 cm Cuvettes	50 PK
MIR 50124	Cell Droppers	25 PK
MIR 50125	Cell Droppers	50 PK

<sup>a</sup> Electroporations per kit.

<sup>b</sup> Number of electroporations in 0.4 cm cuvette.

To inquire about bulk pricing, please call 888.530.0801.

**siRNA****Label IT® RNAi DELIVERY CONTROLS**

- ▷ **Sensitive**—Easily visualize transfected cells and assess delivery efficiency using fluorescent microscopy
- ▷ **Inert**—Does not target any known mammalian genes or cause off-target effects
- ▷ **Compatible**—Suitable for co-delivery experiments with functional siRNA
- ▷ **Easy to Use**—Supplied as a ready to use 10  $\mu$ M stock with a 10X RNAi Dilution Buffer


**PRODUCT SPECIFICATIONS****Label IT® RNAi Delivery Controls**

LABEL	PRODUCT NO.	QUANTITY
Cy™3	MIR 7900	10 $\mu$ g
	MIR 7901	100 $\mu$ g
Fluorescein	MIR 7902	10 $\mu$ g
	MIR 7903	100 $\mu$ g

To inquire about bulk pricing, please call 888.530.0801.

**Description**

The *Label IT* RNAi Delivery Controls consist of either Cy™3 or fluorescein labeled RNA duplex that has the same length, charge and configuration as standard siRNA. The sequence of the *Label IT* RNAi duplex is inert and is not known to affect any cellular events. These controls are designed to facilitate assessment of delivery efficiency of dsRNA oligonucleotides in both *in vitro* and *in vivo* applications and can be co-transfected with a functional target gene-specific siRNA.

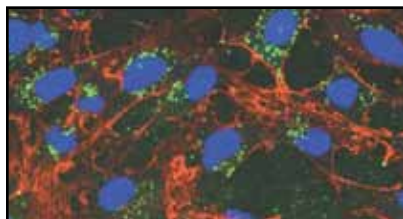

**FIGURE 15.** Visualization of *Label IT*® RNAi Delivery Control. HeLa cells were transfected in complete media with the *Label IT*® Fluorescein RNAi Delivery Control (green) using the *TransIT*®-TKO Transfection Reagent. Twenty-four hours post-transfection, the cells were fixed, then counterstained to locate the nuclei (blue) and actin (red).


## Large RNA (Viral RNA and mRNA)

# TransIT<sup>®</sup>-mRNA TRANSFECTION KIT

- ▷ **High Efficiency RNA Delivery**—Ensures experimental success by effectively transfecting RNA into a large percentage of the cell population
- ▷ **Serum Compatible**—
  - ▶ Simplifies the transfection protocol
  - ▶ Saves valuable time
  - ▶ Increases cell viability thereby minimizing experimental bias due to alterations in cell physiology
- ▷ **Transfects Various RNA Types**—Ideal for multiple applications such as viral production and replication studies and short-term protein expression from mRNA

### Description

*TransIT*-mRNA Transfection Kit provides high efficiency transfections of large RNA molecules such as mRNA or viral RNA. This kit is extremely easy to use and minimizes cellular toxicity due to its unique ability to transfect RNA in the presence of serum.

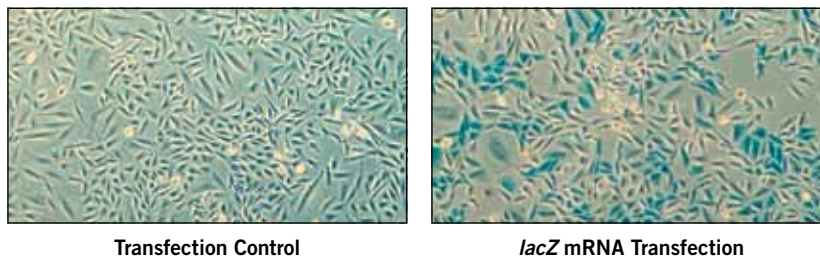


**FIGURE 16.** The *TransIT*<sup>®</sup>-mRNA Transfection Kit Efficiently Delivers *lacZ* mRNA to CHO-K1 Cells. CHO-K1 cells were transfected with an *in vitro* synthesized *lacZ* mRNA using the *TransIT*<sup>®</sup>-mRNA Transfection Kit. Cells were stained 18 hours post-transfection with the Beta-Gal Staining Kit to detect *lacZ* expressing cells.

**FIGURE 17.** The *TransIT*<sup>®</sup>-mRNA Transfection Kit Efficiently Delivers mRNA to a Variety of Cell Lines. Cells were transfected with a capped and polyadenylated EGFP encoding mRNA using the *TransIT*<sup>®</sup>-mRNA Transfection Kit and analyzed by flow cytometry 18 hours post-transfection to identify the EGFP expressing cells.


### PRODUCT SPECIFICATIONS

PRODUCT NO.	QUANTITY
MIR 2225	0.25 ml
MIR 2250	0.5 ml
MIR 2255	5 × 0.5 ml
MIR 2256	10 × 0.5 ml

To inquire about bulk pricing, please call 888.530.0801.

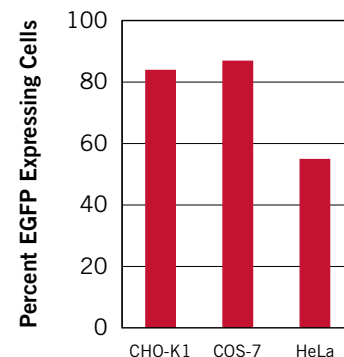

## TransIT<sup>®</sup>-mRNA continued

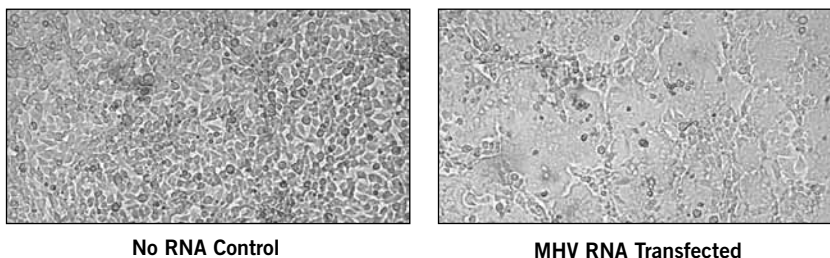

**FIGURE 18.** The *TransIT*<sup>®</sup>-mRNA Transfection Kit Successfully Delivers Viral RNAs at Least 32 kb Long. A 32 kb *in vitro* transcript of the murine coronavirus, MHV, was transfected into DBT cells using the *TransIT*-mRNA Transfection Kit. Successful transfection was assessed by the formation of syncytia 24-48 hours post-transfection. Syncytia were visualized by phase contrast microscopy.


Data courtesy of Mark Clementz, Loyola University of Chicago.

**TABLE 2.** Viral RNAs Successfully Transfected using the *TransIT*<sup>®</sup>-mRNA Transfection Kit.

Virus	Application	RNA Size	Laboratory
Feline Calicivirus	Virus Production	8 kb	Hardy Lab Montana State University
Poliovirus	Virus Production	7.5 kb	Ehrenfeld Lab LID, NIAID, NIH
Hepatitis C Virus	Replicon	9 kb	Striker Lab University of Wisconsin
	Virus Production	9 kb	Striker Lab University of Wisconsin
Yellow Fever Virus	Replicon	6 kb	Striker Lab University of Wisconsin
	Virus Production	6 kb	Striker Lab University of Wisconsin
Murine Coronavirus	Virus Production	32 kb	Baker Lab Loyola University of Chicago

**FIGURE 19.** The *TransIT*<sup>®</sup>-mRNA Transfection Kit Outperforms Electroporation and Saves Valuable Time. 10X Better Performance Using 2.5X Less RNA. Huh7 cells were transfected with 2 µg of a yellow fever virus-luciferase expressing RNA replicon (YFV-luc) using the *TransIT*<sup>®</sup>-mRNA Transfection Kit or electroporated in parallel with 5 µg of the YFV-luc RNA replicon using an ECM630 Electroporator (BTX) in PBS. Twenty-four hours later, cells were harvested and assayed for luciferase activity and total protein content. Luciferase activity was normalized to total protein content to correct for differences in cell number between samples.\*

\* Gonzalez G et al. Selection of an Optimal RNA Transfection Reagent and Comparison to Electroporation for the Delivery of viral RNA. *J Virol Meth* 2007; 145:14.

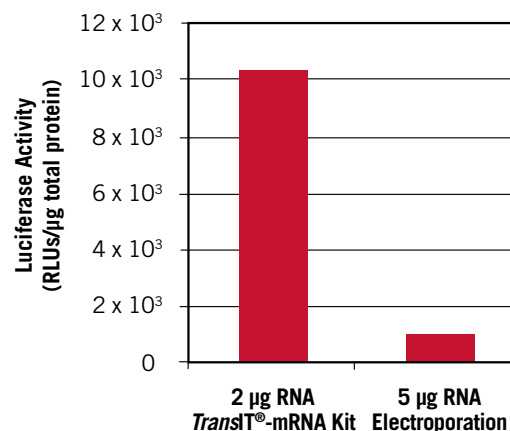





## TERMS AND CONDITIONS

1. **Governing Provisions:** All Mirus Bio LLC products are sold and shipped subject to these Terms and Conditions, unless modified by a quotation issued by Mirus Bio LLC or an agreement signed by both Mirus Bio LLC and the customer. These Terms and Conditions supersede any other written document or oral discussion including terms or conditions appearing on a purchase order issued by the customer.

2. **Product Safety:** Mirus Bio LLC manufactures products for research use by qualified individuals. These products may contain chemicals that may be harmful if misused. Customers are responsible for reviewing and following the instructions included in the product protocols and MSDS that accompany the product. Please contact [techsupport@mirusbio.com](mailto:techsupport@mirusbio.com) with any questions regarding the safe use of Mirus Bio LLC products. We recommend our products be used in accordance with the NIH guidelines developed for genetic research.

3. **Warranties:** Mirus Bio LLC products are warranted to meet or exceed the stated specifications for the usable life of the product defined as either the expiration date for the product or six months from the date of receipt for those products not carrying an expiration date. This warranty is void if the customer has altered or misused the product or failed to store the product as recommended. The warranty does not include defects or failures associated with materials supplied to Mirus Bio LLC by the customer. Mirus Bio LLC disclaims any and all responsibility for any injury or damage which may be caused by the failure of the buyer or any other person to use these products in accordance with the conditions outlined herein. Mirus Bio's sole obligation and the customer's sole remedy are limited to free replacement of products that fail to meet our minimum performance specifications. THIS WARRANTY IS EXCLUSIVE. MIRUS BIO LLC SPECIFICALLY DISCLAIMS ANY OTHER WARRANTIES OF ANY KIND OR NATURE, DIRECT OR INDIRECT, EXPRESS OR IMPLIED, INCLUDING, WITHOUT LIMITATION, THE SUITABILITY, FITNESS OR MERCHANTABILITY FOR ANY PURPOSE. MIRUS BIO WILL NOT BE LIABLE FOR ANY CLAIMS OR DAMAGES UNDER ANY LEGAL THEORY INCLUDING, BUT NOT LIMITED TO CONTRACT, NEGLIGENCE, STRICT LIABILITY OR TORT IN CONNECTION WITH THE FAILURE OF MIRUS BIO LLC PRODUCTS TO PERFORM IN ACCORDANCE WITH THE STATED SPECIFICATIONS. Mirus Bio LLC is not liable for any loss, damages or penalties resulting from delays in manufacturing or delivery of products or services. This warranty applies to all products and services offered by Mirus Bio LLC including catalog products, bulk products, custom products and custom services. Unless otherwise agreed, all technical assistance and information is provided free of charge and we make no warranty regarding the accuracy or utility of such information or assistance.

4. **Patents:** Mirus Bio's products and processes are covered by one or more patents and are subject to other trade secret and proprietary rights. No transfer or grant of rights under any patent or patents is made or is to be implied by any provision under these Terms and Conditions other

than a limited license to use the products for research only. You agree not to infringe upon such rights or to attempt to reverse engineer, reconstruct, synthesize or otherwise modify Mirus Bio products. Certain applications of Mirus Bio products may require licenses from other parties. Determining the existence and scope of such third party intellectual property is the responsibility of the customer. Purchase of the product provides the customer with a limited non-transferable license under any Mirus Bio patents or patent applications to use the product for internal research unless there is a written limitation to this license in the product literature. The customer is responsible for carefully reviewing the product literature and respecting any limitations to this license, e.g. limitations for commercial use or research by for-profit institutions.

5. Customer agrees to indemnify and hold Mirus Bio LLC harmless from and against any and all claims, losses, damages, costs and expenses (including reasonable attorneys' fees and amounts paid in settlement in good faith) which may be suffered or incurred by Mirus Bio LLC (including, without limitation, as a result of any claim or action by a third party) arising in any manner from Customer's use of any Products sold from this Web site or otherwise arising out of any act or omission of Customer, its employees, agents or representatives, or by reason of breach of or failure to perform any of Customer's obligations hereunder.

6. **Orders, Shipment and Delivery:** Purchase orders placed before 4:00 p.m. Central Time will be shipped that day. Orders placed on Friday will ship on the following Monday for delivery on Tuesday. Orders for regularly stocked items may only be cancelled prior to shipment of the order. Shipping charges are prepaid and added to the customer's invoice. Orders are shipped by Federal Express. Unless agreed upon in writing, all shipments are free carrier (FCA) our facility in Madison, Wisconsin. Costs for transportation and risk of loss transfer to the buyer upon delivery to the carrier. The customer is responsible for any tariffs, duties or taxes associated with international shipments.

7. **Copyright and Trademarks:** All content contained in printed materials or on the Mirus Bio LLC website is the copyright of Mirus Bio LLC unless specifically identified otherwise. Trademarks referenced herein are either registered trademarks of Mirus Bio LLC or third party trademarks, which are the property of their respective owners. All rights are reserved. Please contact [sales@mirusbio.com](mailto:sales@mirusbio.com) to request permission to use copyrighted material or trademarks of Mirus Bio LLC.

8. **Payment Terms:** All sales are net 30 days of invoice in U.S. dollars unless other terms are agreed to in writing by Mirus Bio LLC. Prices are subject to change without notice. Mirus Bio LLC accepts the following credit cards: VISA, MasterCard, American Express and Discover. Mail payments to:

**Mirus Bio LLC**  
545 Science Dr.  
Madison, WI 53711 USA  
Attention: Accounts Payable  
Fax: 608.441.2849

9. **Returns:** Mirus Bio LLC must authorize all product returns. Customers wishing to return a product should contact us at [techsupport@mirusbio.com](mailto:techsupport@mirusbio.com). Product that we confirm damaged, incomplete or do not meet our minimum written specifications will be replaced per our warranty (see above) providing such notice is received from the original purchaser within 30 days of receipt of the product. Customers may contact Mirus Bio LLC to cancel an order after shipment or return a product that meets our minimum written specifications. Mirus Bio LLC must authorize these returns. Products must be unopened and returned in their original, intact condition. The customer must bear the cost of shipping the product back to Mirus Bio LLC under the conditions specified by Mirus Bio LLC. All such returns will be subject to a 25% restocking fee.

10. These Terms and Conditions constitute the entire agreement between Mirus Bio LLC and Customer with respect to Customer use of the Web site and Customer purchase of products hereunder, except as the foregoing (i) may be amended from time to time by Mirus Bio LLC, or (ii) as related to Customer purchase of products, may be superseded by any express conflicting terms or supplemented by any express additional terms in a separate written contract signed by authorized representatives of Mirus Bio LLC and Customer.

## PATENT AND LICENSING INFORMATION

Mirus Bio *TransIT*<sup>®</sup> polyamine transfection reagents are covered by U.S. Patent No. 5,744,335, No. 6,180,784, and No. 6,458,382. Mirus Bio *Label IT*<sup>®</sup> nucleic acid labeling and modifying reagents are covered by U.S. Patent No. 6,262,252, and No. 6,593,465. *Cy*<sup>™</sup>3 and *Cy*<sup>™</sup>5 products or portions thereof are manufactured under license from Carnegie Mellon University and are covered by U.S. Patent No. 5,268,486.

Product is sold to the Buyer with a limited license to use the product for research only. The product, or parts from the product, may not be re-packaged or re-sold without written permission from Mirus Bio LLC. A license from Mirus Bio LLC is required for commercial application of the product. For obtaining a license to use the product for commercial application, contact Mirus Bio LLC, 545 Science Drive, Madison, WI 53711. Email: [license@mirusbio.com](mailto:license@mirusbio.com)

## TRADEMARKS

©2010. All rights reserved. Mirus Bio LLC. HeLaMONSTER, HybQUEST, *Label IT*, MiracLEAN,  $\mu$ Array, pLIVE, siQUEST, siXpress, *TransIT-Neural*, *TransIT-TKO* and *TransIT* are registered trademarks of Mirus Bio LLC. Ingenio, Insecta, LOAD 'n GLO and Tracker are trademarks of Mirus Bio LLC.



## DNA TRANSFECTION

### Broad Spectrum

Product Name	Product No.	Quantity
TransIT <sup>®</sup> -2020 Transfection Reagent	MIR 5400	1 ml
	MIR 5404	0.4 ml
	MIR 5405	5 x 1 ml
	MIR 5406	10 x 1 ml
TransIT <sup>®</sup> -LT1 Transfection Reagent	MIR 2300	1 ml
	MIR 2304	0.4 ml
	MIR 2305	5 x 1 ml
	MIR 2306	10 x 1 ml

### Cell Line Specific DNA Transfection

Product Name	Product No.	Quantity
TransIT <sup>®</sup> -293 Transfection Reagent	MIR 2700	1ml
	MIR 2704	0.4 ml
	MIR 2705	5 x 1 ml
	MIR 2706	10 x 1 ml
	MIR 2180	1 ml
TransIT <sup>®</sup> -3T3 Transfection Kit*	MIR 2184	0.4 ml
	MIR 2185	5 x 1 ml
	MIR 2186	10 x 1 ml
	MIR 2170	1 ml
TransIT <sup>®</sup> -CHO Transfection Kit*	MIR 2174	0.4 ml
	MIR 2175	5 x 1 ml
	MIR 2176	10 x 1 ml
	MIR 2190	1 ml
TransIT <sup>®</sup> -COS Transfection Kit*	MIR 2194	0.4 ml
	MIR 2195	5 x 1 ml
	MIR 2196	10 x 1 ml
	MIR 2900	1 ml
TransIT-HeLaMONSTER <sup>®</sup> Transfection Kit*	MIR 2904	0.4 ml
	MIR 2905	5 x 1 ml
	MIR 2906	10 x 1 ml
	MIR 2120	1 ml
	MIR 2124	0.4 ml
TransIT <sup>®</sup> -Jurkat Transfection Reagent	MIR 2125	5 x 1 ml
	MIR 2126	10 x 1 ml
	MIR 2800	1 ml
	MIR 2804	0.4 ml
TransIT <sup>®</sup> -Keratinocyte Transfection Reagent	MIR 2805	5 x 1 ml
	MIR 2806	10 x 1 ml
	MIR 2140	1 ml
	MIR 2144	0.4 ml
TransIT-Neural <sup>®</sup> Transfection Reagent	MIR 2145	5 x 1 ml
	MIR 2146	10 x 1 ml
	MIR 2130	1 ml
	MIR 2134	0.4 ml
TransIT <sup>®</sup> -Prostate Transfection Kit*	MIR 2135	5 x 1 ml
	MIR 2136	10 x 1 ml

### Electroporation

Product Name	Product No.	Size
Ingenio <sup>™</sup> Electroporation Kits (solution, 0.4 cm cuvettes cell droppers)	MIR 50113	25 RXN
	MIR 50116	50 RXN
	MIR 50119	10 RXN
Ingenio <sup>™</sup> Electroporation Kits (solution, 0.2 cm cuvettes cell droppers)	MIR 50112	25 RXN
	MIR 50115	50 RXN
	MIR 50118	10 RXN
Ingenio <sup>™</sup> Electroporation Solution	MIR 50111	25 RXN (6.25 ml)
	MIR 50114	50 RXN (12.5 ml)
	MIR 50117	10 RXN (25 ml)
Ingenio <sup>™</sup> Electroporation Accessories	MIR 50120	0.2 cm cuvettes (25 PK)
	MIR 50121	0.2 cm cuvettes (50 PK)
	MIR 50122	0.4 cm cuvettes (25 PK)
	MIR 50123	0.4 cm cuvettes (50 PK)
	MIR 50124	Cell Droppers (25 PK)
	MIR 50125	Cell Droppers (50 PK)

### NEW! Protein Production

Product Name	Product No.	Quantity
TransIT-PRO <sup>™</sup> Transfection Kit	MIR 5700	1 ml
	MIR 5760	10 ml

### Labeled Plasmid Delivery Controls

Product Name	Product No.	Quantity
Label IT <sup>®</sup> Plasmid	MIR 7904	10 µg
Delivery Control, Cy <sup>™</sup> 3	MIR 7905	100 µg
Label IT <sup>®</sup> Plasmid	MIR 7906	10 µg
Delivery Control, Fluorescein	MIR 7907	100 µg

## RNA TRANSFECTION

### siRNA & miRNA Transfection

Product Name	Product No.	Quantity
TransIT-TKO <sup>®</sup> Transfection Reagent	MIR 2150	1 ml
	MIR 2154	0.4 ml
	MIR 2155	5 x 1 ml
	MIR 2156	10 x 1 ml
	MIR 2110	1 ml
TransIT-siQUEST <sup>®</sup> Reagent	MIR 2114	0.4 ml
	MIR 2115	5 x 1 ml
	MIR 2116	10 x 1 ml
TransIT <sup>®</sup> -siPAK Kit		0.1 ml of each TransIT-siQUEST <sup>®</sup> and TransIT-TKO <sup>®</sup> Reagents
	MIR 2260	
TransIT <sup>®</sup> -siPAK Plus Kit		0.1 ml of each TransIT-siQUEST <sup>®</sup> and TransIT-TKO <sup>®</sup> Reagents and 10 µg Label IT <sup>®</sup> RNAi Delivery
	MIR 2270	Control, Fluorescein

### Electroporation

Product Name	Product No.	Size
Ingenio <sup>™</sup> Electroporation Kits (solution, 0.4 cm cuvettes cell droppers)	MIR 50113	25 RXN
	MIR 50116	50 RXN
	MIR 50119	10 RXN
Ingenio <sup>™</sup> Electroporation Kits (solution, 0.2 cm cuvettes cell droppers)	MIR 50112	25 RXN
	MIR 50115	50 RXN
	MIR 50118	10 RXN
Ingenio <sup>™</sup> Electroporation Solution	MIR 50111	25 RXN (6.25 ml)
	MIR 50114	50 RXN (12.5 ml)
	MIR 50117	10 RXN (25 ml)
Ingenio <sup>™</sup> Electroporation Accessories	MIR 50120	0.2 cm cuvettes (25 PK)
	MIR 50121	0.2 cm cuvettes (50 PK)
	MIR 50122	0.4 cm cuvettes (25 PK)
	MIR 50123	0.4 cm cuvettes (50 PK)
	MIR 50124	Cell Droppers (25 PK)
	MIR 50125	Cell Droppers (50 PK)

### Labeled RNAi Delivery Controls

Product Name	Product No.	Quantity
Label IT <sup>®</sup> RNAi	MIR 7900	10 µg
Delivery Control, Cy <sup>™</sup> 3	MIR 7901	100 µg
Label IT <sup>®</sup> RNAi	MIR 7902	10 µg
Delivery Control, Fluorescein	MIR 7903	100 µg

### mRNA Transfection

Product Name	Product No.	Quantity
TransIT <sup>®</sup> -mRNA Transfection Kit	MIR 2225	0.25 ml
	MIR 2250	0.5 ml
	MIR 2255	5 x 0.5 ml
	MIR 2256	10 x 0.5 ml

MIRUS BIO LLC | 545 Science Dr., Madison, WI 53711  
toll free 888.530.0801 | direct 608.441.2852

Mirus Bio products are distributed worldwide. For a complete list of distributors visit: [www.mirusbio.com/contact/distributors.aspx](http://www.mirusbio.com/contact/distributors.aspx)

Mirus Bio products are also distributed in the U.S. by:



For customer service, call 1-800-766-7000.  
To fax an order, use 1-800-926-1166.  
To order online: [www.fishersci.com](http://www.fishersci.com)